

## Hydrogen Peroxide Breakthrough Indicator Sticker H<sub>2</sub>O<sub>2</sub> BTIS (PN: 1400)

Manual



### 1. Application

The Hydrogen Peroxide BTIS (PN: 1400) is qualitative (yes/no) colorimetric indicator for the saturation and end-of-service life of ductless hood filters. The indicator is designed to provide real-time indication of the breakthrough of hydrogen peroxide.

### 2. Specifications

- |                              |  |
|------------------------------|--|
| a. Weight:                   | 0.4g (0.02oz)                            |
| b. Dimensions:               | 2.8mm (0.11in), $\phi$ : 31.8mm (1.25in) |
| c. Operating temperature:    | -20°C to 45°C (-4°F to 113°F)            |
| d. Operating humidity:       | 5% RH to 85%RH                           |
| e. Minimum detectable limit: | 1ppm-hr at 30 cm/sec face velocity       |
| f. Color change:             | Off white to gray to black.              |
| g. Storage temperature:      | 4°C to 25°C, (39°F to 77°F)              |
| h. Shelf life:               | 1 year at 4°C to 25°C, (39°F to 77°F)    |
| i. Service life:             | 1 year.                                  |

Cross interferences: Strong oxidizers. No other interferences or limitations are known.

### 3. Operating Instructions

- Ensure that packaging pouch is intact.
- Open packaging pouch by tearing off the top part from one of side notches
- Remove indicator sticker from the packaging pouch.
- Peel off the protective liner to expose the bottom adhesive (Figure 1).



**Caution: Do not touch bottom adhesive or exposure area**

- Hold the sticker from the edges, as shown in Figure 2, and place it on center clean area of the filter's outlet with the reading area (glossy surface) of the sticker facing up.
- Press firmly to attach sticker to the filter's outlet (Figure 3).
- place filter when the reading area of the indicator changes color to gray or black.

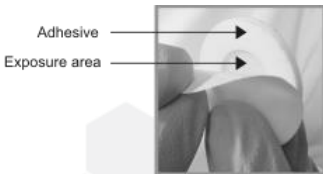


Figure 1



Figure 2

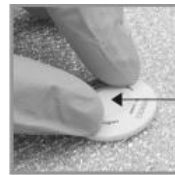


Figure 3

**CHEMTEQ**<sup>®</sup>

Chemteq, Inc., 600 West 24<sup>th</sup> Street, Suite B, Norfolk, Virginia 23517, Tel 757-622-2223, sales@chemteq.net  
©2009 Chemteq, Inc.

[www.chemteq.net](http://www.chemteq.net)

MP-14008100-040424-A